

Installation instruction for Eyespray eye/face wash

Each eyewash station contains of the following items

- 1 pcs rack
- 1 pcs refill with 2 spray bottles,
- 1 pcs red sealing tape and
- 1 pcs white label for notification of the expire date on the bottles.

Art. nr: 822001-3

Mount the rack on a wall.

The rack should be mounted on the height of 1.150 mm above the floor measured up to the bottom of the rack.

So when the eyewash is in an activated position the spray bottles will be on the same height as on a traditional eyewash which is 1.100 mm above the floor.

For installation of the spray bottles see attached: ***Bottle replacement instruction.***

ATTENTION

The red flag indicates that the eyewash has been activated and that new bottles must be installed.

When installing or replacing old spray bottles, ensure that the cover is not completely open so as to avoid inadvertently activating the new bottles.

One way to prevent such accidental activation during the replacement process is to place an object, such as a pencil, between the cover and the backside of the rack.

Bottle replacement instruction

Remove the old spray bottles from the rack. Take two new bottles and note the expire date on the white label. The date is located on the bottom of each bottle.

Remove the cover on each spray bottle.

Install the new bottle with the activator turned towards your chest.

Turn each bottle so the inserter on each bottle is aiming up toward your face.

Simultaneously close the cover and bend the red flag under the cover.

Place the red seal over the edge of the cover and the backside and put, the white label with the expire date on the spray bottles, on the front as attached picture show.



If you need any further assistance please give us a call or send us an e-mail.